

Standard protocols



Protocols

Introduction

Protocols

References

Introduction

These protocols are compiled from the literature from Biacore AB.
These protocols apply to the Biacore 2000/3000 instruments and software

The following protocols are present:


Startup	Starting a Biacore machine
Continue	Leaving an instrument for 3-4 days
Desorb	Cleaning the instrument once a week
Sanitize	Cleaning the instrument once a month
Close	Leaving the instrument up to 5 days
Shutdown	Leaving the instrument for more than 5 days
System check	Checking of the system performance
Emergency stop	Stopping a program at once
Power failure	Recover from power failure

25 February 2016



STARTUP

Start a Biacore 1000, 2000, 3000 instrument. The procedure takes about 1 hour to complete.

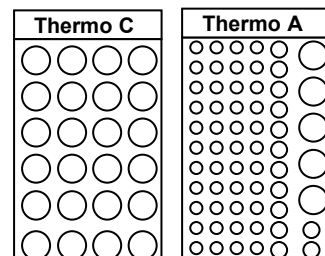

Position of vials

Step / Command	Position	Solution	Volume (ml)	Time (min)	Remarks
Switch on BIAcore instrument	-	-	-	-	-
Wait until the GREEN ready LED light	-	-	-	2	-
Switch on the computer and start BIAcore Control Software	-	-	-	5	-
Make sure there is a Sensor chip inside	-	-	-	-	-
Place buffer in the Bottle compartment and insert tubing	Bottle compartment	Flow buffer	-	-	-
Prime the system	-	Flow buffer	-	7	-
Run Continue/Standby	-	Flow buffer	-	-	max 72 - 96 hours
Wait until the temperature LED is steady	-	-	-	45-60	-

Solution	Preparation
Flow buffer	10 mM HEPES pH 7.4, 150 mM NaCl, 3.4 mM EDTA, 0.005% P20


Standard protocols



Protocols

CONTINUE

Leaving the instrument up to 72 – 96 hours.
 For Control Software 2.x up to 72 hours.
 For Control Software 3.x up to 96 hours.

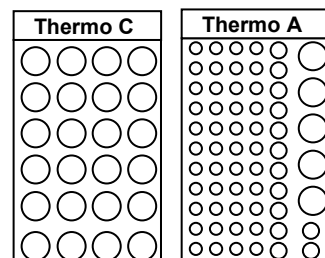

Position of vials


Step / Command	Position	Solution	Volume (ml)	Time (min)	Remarks
Continue/Standby	Bottle compartment	Flow buffer	90	-	After 72 -96 hours the method will stop.

Solution	Preparation
Flow buffer	10 mM HEPES pH 7.4, 150 mM NaCl, 3.4 mM EDTA, 0.005 % P20



DESORB

Clean tubing, IFC-channels and Injection port. Should be done once a week. Use a special desorb chip since the chemicals used can destroy the ligand surface.

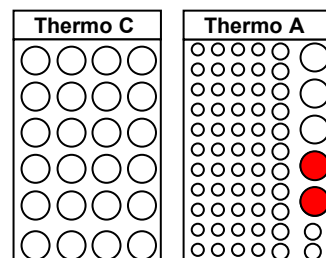

Position of vials

Step / Command	Position	Solution	Volume (ml)	Time (min)	Remarks
Insert desorb chip	-	-	-	-	skip prime
Unclogging	-	Flow buffer	-	4	-
Desorb	R2 F3	Desorb solution 1	3	22	-
	R2 F4	Desorb solution 2	3		
Prime	-	Flow buffer	-	7	-

Solution	Preparation
Desorb solution 1	0.5% (w/v) SDS in water / BIA maintenance kit
Desorb solution 2	50 mM Glycine-NaOH pH 9.5 / BIA maintenance kit
Flow buffer	10 mM HEPES pH 7.4, 150 mM NaCl, 3.4 mM EDTA, 0.005% P20


Standard protocols



Protocols

SANITIZE

Clean tubing, IFC-channels and Injection port from bacteria. Should be done once a month. Use a special desorb chip since the chemicals used can destroy the ligand surface.

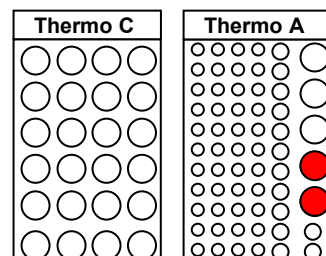

Position of vials

Step / Command	Position	Solution	Volume (ml)	Time (min)	Remarks
Insert Desorb chip	-	-	-	-	skip prime
Flush	-	Flow buffer	-	3	-
Injection port	-	-	-	-	wash injection port by hand with tap water
Needle rep/positioning	-	-	-	-	check needle positions
Unclogging	-	Flow buffer	-	4	-
Desorb	R2 F3 R2 F4	Desorb solution 1 Desorb solution 2	3 3	22	-
Sanitize	Bottle compartment	Hypochlorite sol. Flow buffer	10	35 10	follow messages on the screen
Prime	-	Flow buffer	-	7	-

Solution	Preparation
Desorb solution 1	0.5% (w/v) SDS in water / BIA maintenance kit
Desorb solution 2	50 mM Glycine-NaOH pH 9.5 / BIA maintenance kit
Hypochlorid solution	0.75 ml stock to 10 ml water / BIA maintenance kit
Flow buffer	10 mM HEPES pH 7.4, 150 mM NaCl, 3.4 mM EDTA, 0.005% P20


Standard protocols



Protocols

CLOSE

Leaving the instrument for up to 5 days. Use a special desorb chip since the chemicals used can destroy the ligand surface.

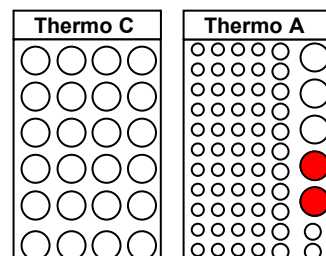

Position of vials

Step / Command	Position	Solution	Volume (ml)	Time (min)	Remarks
Insert Desorb chip	-	-	-	-	skip prime
Unclogging	-	Flow buffer	-	4	-
Desorb	R2 F3	Desorb 1	3	22	-
	R2 F4	Desorb 2	3		
Close	R2F3	water	2.5	11	See messages on screen

Solution	Preparation
Desorb solution 1	0.5% (w/v) SDS in water / BIA maintenance kit
Desorb solution 2	50 mM Glycine-NaOH pH 9.5 / BIA maintenance kit
Flow buffer	10 mM HEPES pH 7.4, 150 mM NaCl, 3.4 mM EDTA, 0.005% P20


Standard protocols

Protocols



SHUTDOWN

Leaving the instrument for more than 5 days. Cleaning tubing, IFC and Injection port, then shutdown the system. Use a special desorb chip since the chemicals used can destroy the ligand surface.
Total time \pm 2 hours.

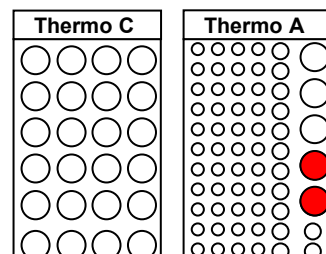

Position of vials


Step / Command	Position	Solution	Volume (ml)	Time (min)	Remarks
Insert Desorb chip	-	-	-	-	skip prime
Change buffer	Bottle compartment	Water	-	3	skip prime
Unclogging	-	-	-	4	-
Desorb	R2 F3 R2 F4	Desorb solution 1 Desorb solution 2	3 3	22	-
Sanitize	Bottle compartment	Hypochlorite sol. Water	10	35 10	follow messages on the screen
Shutdown	Bottle compartment	Water 70 % Ethanol	-	10	follow messages on the screen
Injection port	-	-	-	-	wash injection port by hand with tap water
Needle rep/positioning	-	-	-	-	check needle positions clean vial dislodger
Close the BIACORE Control Software	-	-	-	-	Skip shutdown
Shut off the computer	-	-	-	-	-
Shut off BIACORE	-	-	-	-	-

Solution	Preparation
Desorb solution 1	0.5% (w/v) SDS in water / BIA maintenance kit
Desorb solution 2	50 mM Glycine-NaOH pH 9.5 / BIA maintenance kit
Hypochlorite solution	0.75 ml stock to 10 ml water / BIA maintenance kit
70% ethanol	



SYSTEM CHECK

A comprehensive check of system performance. Insert new CM5 sensor chip first. The procedure must be run at 25°C. Generates a report for evaluation and printing.

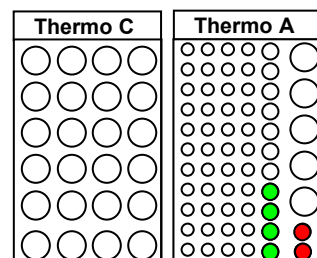

Position of vials


Step / Command	Position	Solution	Volume (ml)	Time (min)	Remarks
Insert new sensor chip	-	-	-	7	Prime
Normalize	R2F2	Glycerol	0.5	5	Run at 25°C
System Check	R2F1	Sucrose	1	40	-
	R2E1	Empty vial	-	-	-
	R2E2	Empty vial	-	-	-
	R2E3	Empty vial	-	-	-
	R2E4	Empty vial	-	-	-
	Bottle Compartment	Flow buffer	-	-	-
System Evaluation	-	-	-	-	-

Solution	Preparation
Glycerol	40% (w/w) Glycerol / BIA maintenance kit
Sucrose	15% (w/w) Sucrose / BIA maintenance kit
Flow buffer	10 mM HEPES pH 7.4, 150 mM NaCl, 3.4 mM EDTA, 0.005% P20



EMERGENCY STOP

Stop a running method at once. Beware that stopping a cycle may result in loss of sample.

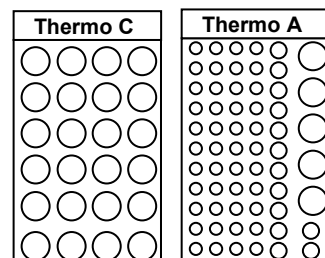

Position of vials


Step / Command	Position	Solution	Volume (ml)	Time (min)	Remarks
Press: Ctrl-BREAK or Run – Stop Method	-	-	-	-	Method will stop when current cycle ends. Read messages.
Press: Ctrl-BREAK again	-	-	-	-	Method will stop immediately. Read messages.
Prime	-	Flow buffer	-	7	To extra clean the system.

Solution	Preparation
Flow buffer	10 mM HEPES pH 7.4, 150 mM NaCl, 3.4 mM EDTA, 0.005% P20



POWER FAILURE

Starting up after a power failure. First try to find the cause !

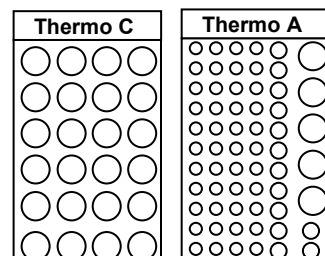

Position of vials

Step / Command	Position	Solution	Volume (ml)	Time (min)	Remarks
First disconnect all plugs or shut down all instruments	-	-	-	-	
Check if the fuse of the power block is on	-	-	-	-	
Then insert BIACORE plug and switch on	-	-	-	-	
Insert the plug of the computer and switch on	-	-	-	-	
Start BIACORE Control Software	-	-	-	-	Save any recovered files first.
When both instruments are working properly, insert the other plugs	-	-	-	7	
Prime	-	Flow buffer	-	-	
Wait until the temperature is stable	-	-	-	-	

Solution	Preparation
Flow buffer	10 mM HEPES pH 7.4, 150 mM NaCl, 3.4 mM EDTA, 0.005% P20

Standard protocols



Protocols

References

References

1. BIACORE AB; BIACORE Application Handbook; 1998
2. BIACORE AB; BIACORE Getting Started; 1998
3. BIACORE AB; BIACORE Instrument Handbook; 1998

Disclaimer

This document is compiled from the publications mentioned above and my own research. The document is provided as is in a good scientific manner to share knowledge.

You may use this document for your own research purposes. You may copy the document as long as you keep it intact. It is not allowed to use this document or parts of it commercially.

For more information, go to the SPR Pages at <http://www.sprpages.nl>.